




VEESPRAY FITTING MAINTENANCE

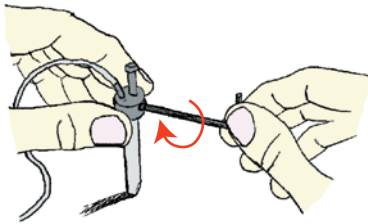


1. Holding the nebulizer in the left hand and leaning it vertically on a smooth flat surface, loosen the 2 grub screws with the right hand, using a small hex key.

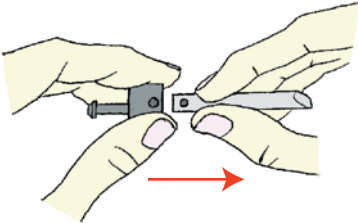

2. Take out the nebulizer body and the old sample tube.

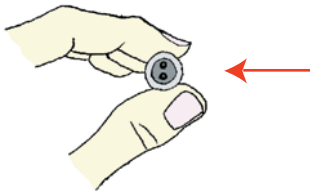


3. Check that the 2 o-rings in the nebulizer fitting are in place.

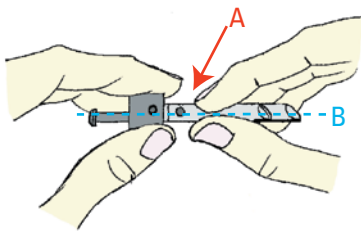


4 a. Replace the nebulizer body into the fitting, lining up the threaded holes with the indentations in the nebulizer body.

b. Line up the gas bore inside the nebulizer body with the gas arm.

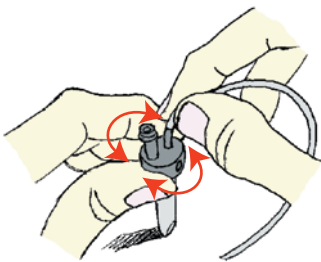

5. Holding the nebulizer in the left hand again, insert the new sample tube by twisting left to right and pushing it in gently, until it stops.

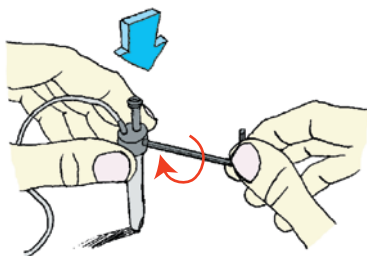

6. Ensure that the indentations in the nebulizer body line up with the threaded holes. Pushing down on the fitting, tighten the 2 grub screws firmly.

Note: Do not use excessive force when tightening.



GLASS EXPANSION

Quality By Design

Head Office (Australia)
15 Batman Street
West Melbourne VIC 3003
Australia
A.B.N. 35 006 560 032

Telephone: +61 3 9320 1111
Toll Free (Aust): 1800 777 638
Facsimile: +61 3 9320 1112
Email: enquiries@geicp.com
www.geicp.com

USA Office
4 Barlows Landing Road
Unit #2
Pocasset, MA 02559
USA

Telephone: 508 563 1800
Toll Free: 800 208 0097
Facsimile: 508 563 1802
Email: geusa@geicp.com